

UN3373 Cat. B

PACKING INSTRUCTION PI 650 IATA

What does the UN3373 kit contain?

Padded bag, absorbent cloth, security bag,
UN3373 Box.



Place the absorbent cloth in the padded bag,
then put your samples in the padded bag as
well. Now put the padded bag in the security
bag and seal the security bag.



Place the security bag in the UN3373 cat B. box
and close the box.




Your sample is now ready for transport.



Goods that are not classified as dangerous goods must be packed in a box or padded envelope.

*If you have any questions about packing, please do not hesitate to contact
CLX - Cargo Logistics Express Tel: 08-594 50062*

DRYICE BOXES

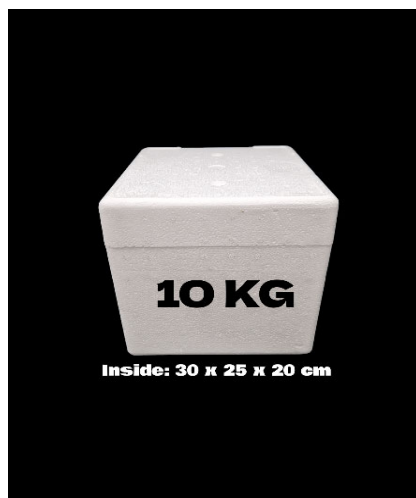

Max Dryice	Size inside
3KG	17x15x14 cm
5KG	25x20x15 cm
10KG	30x25x20 cm

To order dryice and boxes, contact
CLX - Tel: 08-594 50062